

Putting a disinfectant washer out of operation without emptying dispensing system

Instruction:

Putting the dishwasher out of operation without emptying dispensing system

If the machine is not to be used for an extended period of time, please carry-out the following steps.

- Remove all load from machines baskets and inserts
- Remove the filter combination and clean coarse, fine, microfine and flat filter
- Clean the non-return valve and drain pump impeller
- Re-insert the non-return valve and filters
- Clean filters at mobile units accordingly
- Run the Rinsing programme (cold wash cycle) This will ensure that any traces of process are dissolved and rinsed away
- Remove the filter combination and remove residual water using a syringe or sufficient sponge. Finally use a towel for proper drying
- Re-insert the filters
- Close the onsite water intake valves
- Inform the user not to close the machines for at least 12 hours
- Switch the machine off using the On/Off button at the machines control panel
- Switch the machine off at the on-site main switch / fuse breaker

Putting the dishwasher back into operation after a period of non-use

- Open the water intake
- Switch on the on-site main switch/fuse breaker
- Open the machines door
- Check the machine and baskets/mobile units according to the list for daily routine checks
- Check all screw connections in the wash cabinet for proper tightness
- Switch the dishwasher on
- Vent/prime the dispensing system

- Run the Rinsing programme (cold wash cycle) to ensure any traces of process are dissolved
- Run a standard disinfection cleaning programme

